



## Bold SW 1 NV modified medium +2

Roscoff Culture Collection<sup>1</sup>

<sup>1</sup>Station Biologique, CNRS-Sorbonne Université, Roscoff France







ABSTRACT

Medium used for some eukaryotes

Adapted from CCAP web site: https://www.ccap.ac.uk/media/documents/BB.pdf

See also: http://www.ife.ac.uk/ccap/MediaRecipes.html#bb

MATERIALS

NAME CATALOG # **VENDOR** 

View Red Sea Salt Red Sea

BEFORE STARTING

Please refer to our general recommendations to grow cultures :

https://www.protocols.io/private/A48906DC1374AD6281495CB86A8F092F

## Change from standard Bold medium

Volume for 500 ml	Component	Concentration Stock Solutions		
5 ml	CaCl2	CaCl2.2H2O	0.5 gr	200 ml
5 ml	NaNO3	NaNO3	16 gr	200 ml
5 ml	MgSO4	MgSO4.7H2O	10 gr	200 ml
10 ml	P04	From PCRS11 media		
5 ml	NH4Cl (K)	From K media		
3 ml	K trace metal	From K media		
50 μΙ	Vitamines			
467 ml	SW sterile			