

Preparation BG11 medium

1. Under hood, to 1L of sterile water, add 20mL of Cyanobacteria BG-11 Freshwater Solution (Sigma-Aldrich C3061)
2. Filter the medium on 0,2microns

The RCC team



Station Biologique de Roscoff,
CNRS/Université P. & M. Curie,
Place G. Teissier,
29680 Roscoff, France

*Apparatus : Autoclave, Laminar
flow cabinet*

*Solutions : Sterile water,
Nutriments (see protocol)*

*Plasticware and filters : Bottle in
polycarbonate Nalgene, Pipette,
Stericup® Filter Unit (Millipore,
SCGVU10RE)*

ASSEMBLE

ASSOCIATION OF EUROPEAN MARINE BIOLOGICAL LABORATORIES