

Preparation K/2 medium

1. Under hood, to 500mL of seawater beforehand filter and autoclaved, add 500mL of K medium (cf. protocol K medium)
2. Filter the medium on 0,2microns (optional)

The RCC team



Station Biologique de Roscoff,
CNRS/Université P. & M. Curie,
Place G. Teissier,
29680 Roscoff, France

*Apparatus : Autoclave, Laminar
flow cabinet, Stainless Steel
Filter Holders*

Solutions : Seawater, Medium K

*Plasticware and filters : Bottle in
polycarbonate Nalgene, Glass
fibre prefilters (Millipore,
AP1507500), Filters 0,22µm
GSWP (Millipore, GSWP09000),
Stericup® Filter Unit (Millipore,
SCGVU10RE)*

ASSEMBLE

ASSOCIATION OF EUROPEAN MARINE BIOLOGICAL LABORATORIES