

ASW+NO3 medium Roscoff Culture Collection¹¹CNRS-Sorbonne Université, Station Biologique, Place G. Tessier 29680 Roscoff FRANCE

Daniel Vaultot

Station Biologique, Roscoff, France



ABSTRACT

Medium used for cyanobacteria based on artificial seawater (ASW)

BEFORE STARTING

Please refer to our general recommendations to grow cultures :

<https://www.protocols.io/private/A48906DC1374AD6281495CB86A8F092F>

- 1
 - Dissolve 25 g of NaCl in MilliQ water
 - To this solution, add :

<i>Quantity</i>	<i>Compound</i>	<i>Stock Solution</i>	<i>Concentration in medium (in mM)</i>	
<i>mLstock /LASW</i>	<i>g/LASW</i>			
10	0,75	Sodium nitrate (NaNO ₃)	75 g/L	8.8
10	2	Magnesium chloride hexahydrate (MgCl ₂ 6H ₂ O)	200 g/L	9.8
5	0,5	Potassium chloride (KCl)	100 g/L	6,7
5	0,5	Calcium chloride (CaCl ₂)	100 g/L	4.5
10	3,5	Magnesium sulfate heptahydrate (MgSO ₄ 7H ₂ O)	350 g/L	14.2
5,5	1,1	TRIS-Base	200 g/L	9.08
2,5	0,03	Dipotassium phosphate (K ₂ HPO ₄)	12 g/L	0.172

- Adjust the pH to 8 with concentrated HCl
- Adjust to 999 mL with milliQ water
- Add 1 mL of trace metals (see receipe below)
- Autoclave the medium

Trace metal stock solution

- 2
- Dissolve all these components separately in milliQ water :

<i>Quantity</i>	<i>Compound</i>
2.86g	Boric acid (H ₃ B ₃ O ₃)
1.81g	Manganese (II) chloride tetrahydrate (MnCl ₂ ·4H ₂ O)
0.222g	Zinc sulfate monohydrate (ZnSO ₄ ·H ₂ O)
0.390g	Sodium molybdate dihydrate (Na ₂ MoO ₄ ·2H ₂ O)
0.008g	Copper sulfate pentahydrate (CuSO ₄ ·5H ₂ O)
0.0494g	Cobalt nitrate hexahydrate (Co(NO ₃) ₂ ·6H ₂ O)
3.0g	Ferric chloride hexahydrate (FeCl ₃ ·6H ₂ O)
0.5g	EDTA magnesium disodium (EDTA(Na ₂ Mg))

- Combine the various solutions after full dissolution
- Make final volume up to 1L with milliQ
- Store in refrigerator